Bond™ Intense R Detection

Catalog No: DS9263

Intended Use

For research use only. Not for use in diagnostic procedures.

Bond Intense R Detection is a biotin/streptavidin system for the detection of tissue-bound primary antibodies. It is intended for staining sections of formalin-fixed, paraffin-embedded tissue on the Bond automated system. The user must supply a biotinylated secondary antibody of choice.

Summary and Explanation

Immunohistochemical techniques can be used to demonstrate the presence of antigens in tissue and cells (see "Using Bond Reagents" in your Bond user documentation).

Bond Intense R Detection works as follows:

- The specimen is incubated with hydrogen peroxide to quench endogenous peroxidase activity.
- A user-supplied specific primary antibody is applied.
- A user-supplied biotinylated secondary antibody localizes the primary antibody.
- A streptavidin-enzyme conjugate binds to the biotin present on the secondary antibody.
- The substrate chromogen, 3,3'-diaminobenzidine (DAB), visualizes the complex via a brown precipitate.
- Hematoxylin (blue) counterstaining allows the detection of cell nuclei.

Reagents Provided

1. Peroxide Block (30 mL) 3-4% (v/v) Hydrogen peroxide.
2. Open Secondary Antibody – empty container in which to place the biotinylated secondary antibody of choice.
3. Streptavidin HRP (30 mL) (1-1.4 µg/mL) in Tris buffered saline with protein stabilizer and 0.35% ProClin™ 950.
4. DAB Part 1 (2.4 mL) 66 mM 3,3'-diaminobenzidine tetrahydrochloride hydrate, in a stabilizer solution.
5. DAB Part 2 (30 mL) 9% (v/v) N,N-Dimethylformamide in a stabilizer solution.
6. DAB Part 2 (30 mL) 9% (v/v) N,N-Dimethylformamide in a stabilizer solution.
7. DAB Part 3 (2.4 mL) 0.75% (v/v) Hydrogen Peroxide.
8. Hematoxylin (30 mL) <0.1% Hematoxylin.

Reconstitution, Mixing, Dilution, Titration

Bond Intense R Detection is optimized for use on the Bond system. Reconstitution, mixing, dilution, or titration of these reagents is not required.

Storage and Stability

Store at 2-8 °C. Do not freeze. Do not use after the expiration date indicated on the tray handle label. Return to 2-8 °C immediately after use. Any storage conditions other than those specified above must be verified by the user¹.

Instructions for Use

Fill the container labelled "Open Secondary Antibody" with 35 mL of a biotinylated secondary antibody of choice, diluted in Bond Primary Antibody Diluent.

The optimal concentration and stability of this reagent must be determined by the user.

Bond Intense R Detection was developed for use on the Bond automated system.

Operating parameters for application of the detection system reagents on the Bond Processing Module have been optimized at Leica Biosystems. These can be displayed by following the instructions in your Bond user documentation.

Troubleshooting

Refer to reference 4 for remedial action. Contact your local distributor or the regional office of Leica Biosystems to report unusual staining.

Further Information

Further information on immunostaining with Bond reagents, under the headings Principle of the Procedure, Materials Required, Specimen Preparation, Quality Control, Assay Verification, Interpretation of Staining, Key to Symbols on Labels, and General Limitations can be found in "Using Bond Reagents" in your Bond user documentation.

Bibliography

1. Clinical Laboratory Improvement Amendments of 1988. Final Rule 57 FR 7163 February 28, 1992.
2. List of substances which may be candidates for further scientific review and possible identification, classification and regulation as potential occupational carcinogens. Fed Reg 1980;45:157.
3. Villanova PA. National Committee for Clinical Laboratory Standards (NCCLS). Protection of laboratory workers from infectious diseases transmitted by blood and tissue; proposed guideline. 1991; 7(9). Order code M29-P.
4. Bancroft JD and Stevens A. Theory and Practice of Histological Techniques. 4th Edition. Churchill Livingstone, New York. 1996.

Warnings and Precautions

- To obtain a copy of the Material Safety Data Sheet contact your local distributor or regional office of Leica Biosystems, or alternatively, visit the Leica Biosystems Web site, www.LeicaBiosystems.com
- Specimens, before and after fixation, and all materials exposed to them, should be handled as if capable of transmitting infection and disposed of with proper precautions². Never pipette reagents by mouth and avoid contacting the skin and mucous membranes with reagents or specimens. If reagents or specimens come in contact with sensitive areas, wash with copious amounts of water. Seek medical advice.
- Consult Federal, State or local regulations for disposal of any potentially toxic components.
- Minimize microbial contamination of reagents or an increase in non-specific staining may occur.
- Incubation times or temperatures other than those specified may give erroneous results. Any such change must be validated by the user¹.
- Do not mix reagents from different detection systems.

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© Leica Biosystems Newcastle Ltd · DS9263 02/03/2015
Leica Biosystems Newcastle Ltd