


## **Kreatech™ FISH probes**

### **Product Information Sheet**

KI-60005  
FISH Reagent Kit  
5x5 slides

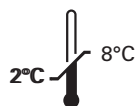



**Kreatech Biotechnology B.V.**  
Vlierweg 20  
1032 LG Amsterdam  
The Netherlands  
[www.LeicaBiosystems.com](http://www.LeicaBiosystems.com)

**RUO - Research Use Only**

Not for use in diagnostic procedures

## Kreatech™ FISH Reagent Kit

### Kit Content

Code #	Description	Volume	Signal Word	Pictogram
LK-095B	DAPI/Antifade 0.1 µg/ml (DAPI Counterstain)	0.5 ml	Warning	
LK-097B	Antifade (Counterstain diluent)	0.5 ml	Warning	
LK-102B	0.4 x SSC / 0.3% Igepal (Wash Buffer I)	250 ml	N.A.	N.A.
LK-103B	2 x SSC / 0.1% Igepal (Wash Buffer II)	2 x 250 ml	N.A.	N.A.
LK-105B	2 x SSC / 0.5% Igepal	250 ml	N.A.	N.A.

### Notes:

- All reagents are ready to use.
- For more intense counterstain DAPI Counterstain 1 µg/mL (LK-096A) is available.

**Warning and precautions:** In case of emergencies check SDS sheets for medical advice. SDS sheets may be obtained by either contacting Leica Technical Support or visiting [www.LeicaBiosystems.com](http://www.LeicaBiosystems.com). DNA probes contain formamide which is a teratogen; do not inhale or allow skin contact. Wear gloves and a lab coat when handling DNA probes. All materials should be disposed of according to your institution's guidelines for hospital waste disposal.

### Reagent Storage and Handling:

Store at 2-8 °C. Reagents should not be used after the expiration date on the vial label.

### TECHNICAL SUPPORT

Technical support is available at [www.LeicaBiosystems.com/service-support/technical-support/](http://www.LeicaBiosystems.com/service-support/technical-support/) or toll free at 800-248-0123 or via e-mail: [kreatech-support@leicabiosystems.com](mailto:kreatech-support@leicabiosystems.com).

### CUSTOMER SERVICE

Kreatech probes may be ordered through Leica Customer Service toll free at 800-248-0123 or order via e-mail: [purchase.orders@leica-microsystems.com](mailto:purchase.orders@leica-microsystems.com).