# Novocastra™ Fluorescein-Conjugated Epstein-Barr virus Probe

BIO SYSTEMS

## Product Code: NCL-EBV

**Intended Use** FOR RESEARCH USE ONLY.

**Product Preparation** Fluorescein labeled oligonucleotide cocktail for the detection of mRNA sequences contained in 1 mL of hybridization solution sufficient for in situ hybridization staining of 50 preparations.

**Storage Conditions** Store at 2–8 °C in the dark. The product is stable under these conditions up to the expiry date indicated on the vial label.

**General Overview** The Epstein-Barr virus (EBV) probe demonstrates cells latently infected with EBV. The probe hybridizes to abundantly expressed Epstein-Barr virus-encoded RNA (EBER) transcripts which are concentrated in the nuclei of latently infected cells. These transcripts are thought to block the activation of dsRNA-dependent eukaryotic initiation factor 2a (eIF-2a) protein kinase DAI. In the absence of EBER, eIF-2a inhibits cellular protein synthesis. The EBER molecules therefore, promote both host cell survival and viral growth (Clarke, Glickman). In situ hybridization using the EBV probe can be undertaken within the working day and completed after overnight substrate incubation. The method is reproducible and results in intense blue/black nuclear staining of cells containing EBV against a virtually clear background. Furthermore, due to the destruction of RNases by formalin fixation, the EBV sequences are conserved in routine laboratory preparations (Pringle).

**General References** Cinque P, Brytting M, et al. *Lancet*. 342: 398–401 (1993).  
Prange E, Trautmann J C, Kreipe H, et al.. *Journal of Pathology*. 166: 113–119 (1992).  
Pringle J H, Barker S, Warford A. *Journal of Pathology (Suppl)*. 167: 133A (1992).  
Sandvej K, Krenacs L, Hamilton-Dutoit S J, et al.. *Histopathology*. 20: 387–395 (1992).  
Wu T-C, Mann R B, Epstein J I, et al.. *American Journal of Pathology*. 138: 1461–1468 (1991).  
Clarke P A, Sharp N A and Clemens M J. *European Journal of Biochemistry*. 193: 635–641 (1990).  
Wu T-C, Mann R B, Charache P, et al.. *International Journal of Cancer*. 46: 801–804 (1990).  
Glickman J N, Howe J G and Steitz J A. *Journal of Virology*. 62: 902–911 (1988).  
Okano M, Thiele G M, Davis J R, et al.. *Clinical Microbiology Reviews*. 1: 300–312 (1988).

## Warnings and Precautions

### EPSTEIN-BARR VIRUS PROBE

Contains Formamide (<30%) and Dextran Sulphate (<10%).  
GHS07: Exclamation mark.

GHS08: Health hazard.

Signal words:  
Danger.

H315: Causes skin irritation.

H319: Causes serious eye irritation.

H360D: May damage the unborn child.

P201: Obtain special instructions before use.

P202: Do not handle until all safety precautions have been read and understood.  
P260: Do not breathe dust/fumes/gas/mist/vapours/spray.

P281: Use personal protective equipment as required.

P308 + P313: IF exposed or concerned: Get medical advice/attention.

P314: Get medical advice/attention if you feel unwell.

P264: Wash thoroughly after handling.

P280: Wear protective gloves/protective clothing/eye protection/face protection.

P305 + P351 + P338: IF IN EYES: Rinse cautiously with water for several minutes.

Remove contact lenses, if present and easy to do. Continue rinsing.

P337 + P313: If eye irritation persists: Get medical advice/attention.

P302 + P352: IF ON SKIN: Wash with plenty of soap and water.

P332 + P313: If skin irritation occurs: Get medical advice/attention.

P362: Take off contaminated clothing and wash before reuse.

Restricted to professional users.

U.S. Patent No. 5,597,692 licensed from Syngene Ltd.

Leica Biosystems Newcastle Ltd  
Balliol Business Park West  
Benton Lane  
Newcastle Upon Tyne NE12 8EW  
United Kingdom  
J +44 191 215 4242

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# Instructions for Use

## In Situ Hybridization Method

Please read the procedural notes before commencing this protocol.

### Sample Preparation

1. Dewax sections in xylene for 2 x 3 minutes.
2. Hydrate in 99% v/v ethanol for 2 x 3 minutes.
3. Hydrate in 95% v/v ethanol for 3 minutes.
4. Immerse in water for 3 minutes (see Procedural Notes, point 4).
5. Place slides on an incubation tray and cover with 100  $\mu$ L of Proteinase K in 0.05 M Tris/HCl buffer pH 7.6 and incubate for 30 minutes at 37 °C.
6. Immerse in water for 2 x 3 minutes (see Procedural Notes, point 4).
7. Dehydrate in 95% v/v ethanol for 3 minutes.
8. Dehydrate in 99% v/v ethanol for 3 minutes.
9. Air dry.

### Hybridization

1. Add 20  $\mu$ L of probe hybridization solution to slides as required and coverslip sections.
2. Incubate for 2 hours at 37 °C (see Procedural Notes, point 6).
3. Allow coverslips to drain off into a beaker.
4. Wash slides in TBS containing 0.1% v/v Triton X-100 for 3 x 3 minutes.

### Detection

Reagents supplied in the Novocastra™ In Situ Hybridization Detection Kit (NCL-ISH-D) can be used in steps 2 and 5 of the following method. When the kit is used, please refer to the supplied instructions for use of the reagents.

1. Place slides on incubation tray and cover sections with 100  $\mu$ L of normal rabbit serum diluted 1:5 in TBS containing 3% w/v BSA, 0.1% v/v Triton X-100. Incubate for 10 minutes.
2. Tip off the blocking solution and add Rabbit F(ab') anti-FITC conjugated to alkaline phosphatase diluted 1:200 in TBS containing 3% w/v BSA and 0.1% v/v Triton X-100. Incubate for 30 minutes.
3. Wash slides in TBS for 2 x 3 minutes.
4. Wash slides in alkaline phosphatase substrate buffer pH 9.0 for 5 minutes.
5. Place slides in incubation tray and demonstrate alkaline phosphatase activity by covering the sections with 100  $\mu$ L of the following solution: 10  $\mu$ L Alkaline Phosphatase substrate buffer, 80  $\mu$ L 5-bromo-4-chloro-3-indolyl phosphate (BCIP) (50 mg/mL in dimethyl formamide), 80  $\mu$ L Nitro blue tetrazolium (NBT) (75 mg/mL in 70% v/v dimethyl formamide), 10  $\mu$ L 1 M levamisole, coverslip and incubate at room temperature in the dark overnight.
6. Wash in running water for 5 minutes.
7. Counterstain in Mayer's hematoxylin for a maximum of 10 seconds.
8. Mount in aqueous mountant.

### Suggested Solutions

**TBS:** 50 mM Tris/HCl, 150 mM NaCl pH 7.6

**Alkaline Phosphatase substrate buffer:** 100 mM Tris/HCl, 50 mM MgCl<sub>2</sub>, 100 mM NaCl pH 9.0.

### Procedural Notes

1. Use of the control probe (NCL-CONTROL) is recommended alongside the application of specific probes to control for non-specific probe interactions with tissue. In the small and large intestine, cells of the diffuse endocrine system will react non-specifically with the probes and in this situation the control probe should always be applied to duplicate sections.
2. For surgical tissues fixed in formalin for 24–48 hours and paraffin wax-embedded, digestion with a range of Proteinase K concentrations of 5, 10 and 15  $\mu$ g/mL should be carried out. Under and over digestion will result in suboptimal signal strength.
3. The probe hybridization solution contains formamide and care should be taken to avoid skin contact.
4. All solutions should be prepared using high grade chemicals and highest quality water available in the laboratory, ie, reverse osmosis and deionized or double distilled.
5. For tissues fixed in solutions containing mercury, it is necessary to pretreat sections before the Proteinase K step with 0.2 M HCl for 20 minutes, then wash in water for 2 x 5 minutes. This procedure is essential to reduce background staining resulting from mercury fixation.
6. Endogenous Alkaline Phosphatase activity can be blocked by denaturing sections prior to hybridization. For denaturation, the preparations should be covered with the probe solution and coverslipped as for the standard procedure. The preparation should then be heated at 65 °C for 15 minutes, after which hybridization at 37 °C should proceed for the standard time of 2 hours.