Novocastra™ Fluorescein-conjugated probes for in situ hybridization ISH Detection Kit

Product Code: NCL-ISH-D

Intended Use

FOR RESEARCH USE ONLY.

Contents

1. Introduction
2. Kit Contents
3. Materials Required but not Provided
4. Preparation of Reagents
5. Precautions and Safety
6. Storage
7. Procedure

Introduction

The Novocastra ISH Detection Kit is intended for the detection of fluorescein-conjugated nucleic acid probes hybridized to their targets in cell or tissue preparations.

The kit can be used in conjunction with the Novocastra range of fluorescein-conjugated probes. Sufficient reagents for the in situ hybridization staining of 100 preparations are provided.

The hybridization is followed by a wash with Tris-buffered saline. Alkaline phosphatase-conjugated Rabbit F(ab') Anti-FITC is then added, and the specimen is incubated at room temperature for 30 minutes followed by a wash. Non-specific binding to the specimen is kept at a minimum due to the use of an antibody F(ab') fragment.

Enzyme substrate (BCIP/NBT) is added, and the specimen incubated at room temperature in the dark overnight (Note after hybridization with the Novocastra Poly d(T) probe, substrate incubation time need not exceed 60 minutes).

Finally the specimen is mounted. Positive staining is recognized under the microscope as a dark blue/black color at the site of hybridization.

Kit Contents

Vial A Rabbit F(ab') Anti-FITC/AP: 0.25 mL

Alkaline phosphatase-conjugated antibody to fluorescein isothiocyanate (affinity-isolated rabbit F(ab')). In 100 mM Tris buffer, 50 mM MgCl₂, 0.1 mM ZnCl₂, 4% w/v BSA, stabilizer, preservative, pH 7.5.

Vial B Enzyme Substrate: 0.4 mL

(50 x concentrated) 5-Bromo-4-Chloro-3-Indolyl Hydrogen Phosphate-P-Toluidine (BCIP) and Nitro Blue Tetrazolium (NBT, Ditetrazolium Dichloride) in Dimethyl Formamide solution.

Vial C Inhibitor (Levamisole): 0.15 mL

(1000x concentrated) 1 M Levamisole Hydrochloride.

Materials Required but not Provided

1. General laboratory equipment for histological procedures.
2. Equipment and reagents needed for the hybridization step.
3. Silanized slides.
4. Tris-buffered saline (TBS): 50 mM Tris/HCl, 150 mM NaCl, pH 7.6.
5. TBS, 0.1% v/v Triton X-100.

6. TBS, 3% w/v BSA, 0.1% v/v Triton X-100, (conjugate diluent).
7. TBS, 3% w/v BSA, 0.1% v/v Triton X-100, 20% v/v normal rabbit serum.
8. Rabbit serum (normal).
9. 100 mM Tris/HCl, 50 mM MgCl₂, 100 mM NaCl pH 9.0 (alkaline phosphatase substrate buffer).
10. Aqueous mountant.

Preparation of Reagents

Anti-FITC/AP Working Solution

Dilute Rabbit F(ab') Anti-FITC/AP (Vial A) 1:100 to 1:200 in TBS, 0.1% v/v Triton X-100, 3% w/v BSA.

Enzyme Substrate with Inhibitor

Dilute Enzyme Substrate (Vial B) 1:50 in 100 mM Tris/HCl, 50 mM MgCl₂, 100 mM NaCl (TBS) pH 9.0. Add 1 µL of Inhibitor (Levamisole) (Vial C) to each mL of diluted Enzyme Substrate.

Storage

The reagents of the kit should be stored in the dark at 2–8 °C.

Procedure

To ensure that sections adhere to the slides throughout the procedure, it is recommended to use silanized slides.

Refer to the in situ hybridization procedure supplied with Novocastra oligonucleotide probes for pretreatment and hybridization steps of the protocol.

Detection

1. Place slides in an incubation tray and cover sections with 100 µL of TBS, containing 3% w/v BSA, 0.1% w/v Triton-X-100 and 20% w/v normal rabbit serum. Incubate for 10 minutes.
2. Tip off the blocking solution and add Rabbit F(ab') Anti FITC/AP diluted 1:100 to 1:200 in TBS, 3% w/v BSA, 0.1% w/v Triton X-100. Incubate for 30 minutes.
3. Wash slides in TBS for 2 x 3 minutes.
4. Wash slides in alkaline phosphatase substrate buffer pH 9.0 for 5 minutes.
5. Place slides in incubation tray and demonstrate alkaline phosphatase activity by covering sections with the following solution: Dilute Enzyme Substrate 1:50 in 100 mM Tris/HCl, 50mM MgCl₂, 100 mM NaCl, pH 9.0. Add 1 µL of Inhibitor (Levamisole) to each mL of diluted Enzyme Substrate. Incubate at room temperature in the dark overnight.
6. Wash in running water for 5 minutes.
7. Counterstain in Mayer's hematoxylin for 10 seconds maximum.
8. Mount in aqueous mountant.

Precautions and Safety

For additional information on toxicity please refer to the Safety Data Sheets (MSDS) for the above components, which are available on request.

Due to the small volumes in the stock reagent sarstedt vials, please ensure that each reagent is in the vial and not the vial cap before opening. This can be achieved simply, by either centrifugation or gently tapping the side of the vial.

Enzyme Substrate

Contains Dimethyl Formamide (<100%), Ditetrazolium Dichloride (<5%) and 5-Bromo-4-Chloro-3-Indolyl Hydrogen Phosphate-P-Toluidine (<5%).
GHS07: Exclamation mark.
GHS08: Health hazard.
Signal words: Danger.

H360D: May damage the unborn child.
H315: Causes skin irritation.
H319: Causes serious eye irritation.
H302+312+332: Harmful if swallowed, in contact with skin or if inhaled.

P201: Obtain special instructions before use.
P202: Do not handle until all safety precautions have been read and understood.
P261: Avoid breathing dust/fumes/gas/mist/vapours/spray.
P263: Avoid contact during pregnancy/while nursing.
P264: Wash hands thoroughly after handling.
P270: Do not eat, drink or smoke when using this product.
P271: Use only outdoors or in a well-ventilated area.
P280: Wear protective gloves/protective clothing/eye protection/face protection.
P301+P312: IF SWALLOWED: Call a POISON CENTRE or doctor if you feel unwell.
P330: Rinse mouth.
P302+P352: IF ON SKIN: Wash with plenty of soap and water.
P332+P313: If skin irritation occurs: Get medical advice/attention.
P304+P340: IF INHALED: Remove victim to fresh air and keep at rest in a position comfortable for breathing.
P305+P351+P338: IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.
P337+P313: If eye irritation persists: Get medical advice/attention.
P308+P313: IF exposed or concerned: Get medical advice/attention.
P312: Call a POISON CENTRE or doctor if you feel unwell.
P362 + P364: Take off contaminated clothing and wash it before reuse.
P501: Dispose of contents/container to hazardous or special waste collection point.
Restricted to professional users.

Inhibitor

Contains Levamisole Hydrochloride (<30%).
GHS07: Exclamation mark.
Signal words: Warning.

H302: Harmful if swallowed.

P264: Wash hands thoroughly after handling.
P270: Do not eat, drink or smoke when using this product.
P301+P312: IF SWALLOWED: Call a POISON CENTRE/doctor if you feel unwell.
P501: Dispose of contents/container to hazardous or special waste collection point.

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Leica Biosystems Newcastle Ltd
Leica Biosystems Inc
Pty Ltd